

## Exchanging the handle of C6100-2G Syringe

1. Pull the plunger out from the syringe




2. Take the end of the piston with the handle as shown here:



Use the index finger and middle finger and hold the white ring on the bottom side - be careful to touch only the white ring, not the black handle.

Now press out the black handle with a little force from your thumb into the indicated direction, the handle should come out.

3. If it is not possible to move the handle through exertion of force another option would be to loosen the handle through knocking carefully on a hard surface.

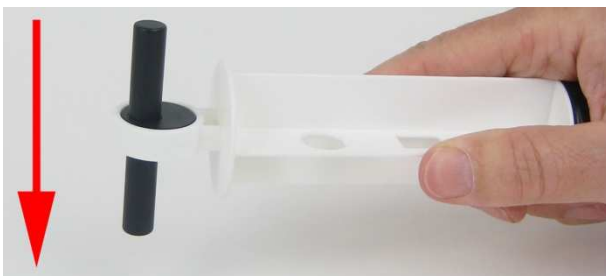

4. To insert the handle into the new syringe, repeat the steps in reverse order.