



**einstein™**  
imagine • explore • learn

## einstein™ Splitter


Product Number: ENSPL011



### Overview

The Einstein Splitter allows you to add even more sensors to your einstein™ Labmate™ or einstein™ Tablet.

### Experiment set up

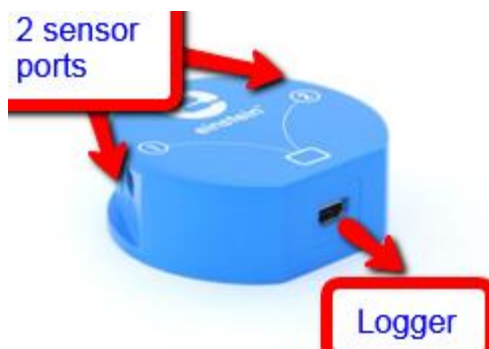

The Splitter fits into one of the mini USB 8 ports located on the side of the data logger. After inserting the splitter two devices can now be plugged into the port.

Note: Each Splitter can support two analog devices or one analog and one digital device but not two digital devices.

## Troubleshooting

If the devices connected to the Splitter aren't automatically recognized by MultiLab/ MiLAB, please contact Fourier Education's technical support.

## Technical support

For technical support, you can contact the Fourier Education's technical support team at:

Web: [www.einsteinworld.com/support](http://www.einsteinworld.com/support)

Email: [support@fourieredu.com](mailto:support@fourieredu.com)

Phone (in the US): (877) 266-4066

## Copyright and Warranty

All standard Fourier Systems sensors carry a one (1) year warranty, which states that for a period of twelve months after the date of delivery to you, it will be substantially free from significant defects in materials and workmanship.

This warranty does not cover breakage of the product caused by misuse or abuse.

This warranty does not cover Fourier Systems consumables such as electrodes, batteries, EKG stickers, cuvettes and storage solutions or buffers.