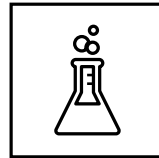
- DE** Bedienungsanleitung
- EN** Operating Instructions
- FR** Mode d'emploi
- NL** Handleiding
- IT** Istruzioni per l'uso
- ES** Instrucciones de uso
- PT** Manual de utilização

MANUAL DOWNLOAD:

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EXPERIMENTS:

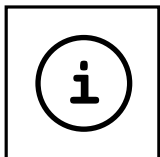
www.bresser.de/downloads

SERVICE AND WARRANTY:

www.bresser.de/warranty_terms

MICROSCOPE GUIDE:

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

MICROSCOPE FAQ:

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DE	Bedienungsanleitung	4
EN	Operating Instructions	12
FR	Mode d'emploi	20
NL	Handleiding	28
IT	Istruzioni per l'uso	36
ES	Instrucciones de uso	44
PT	Manual de utilização	52

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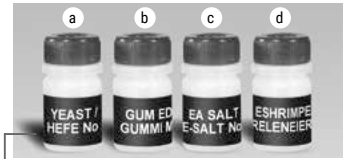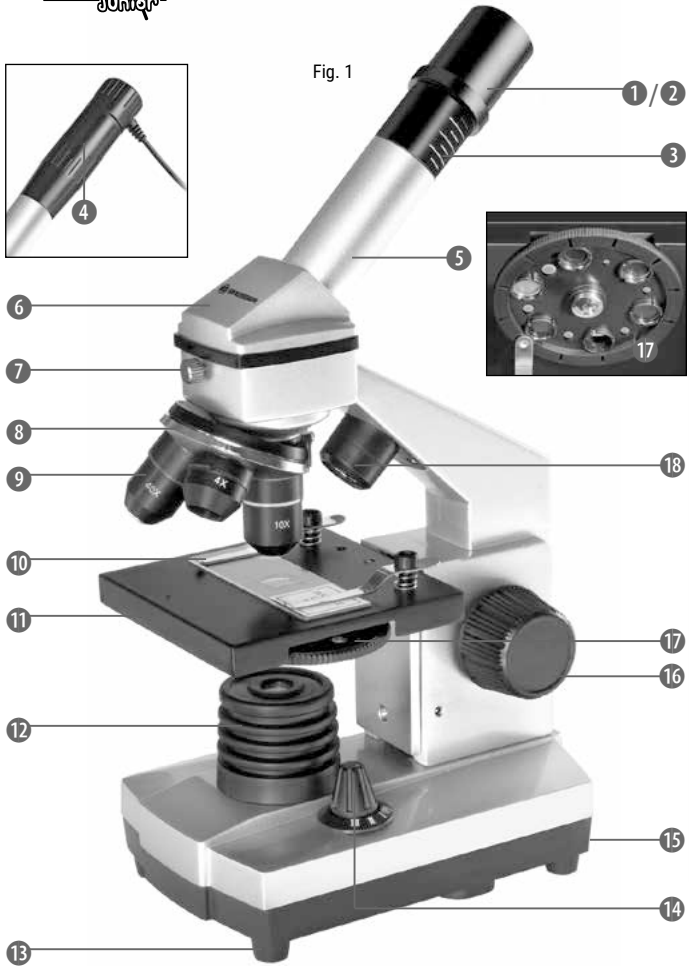RISK to your child.

 Aids with sharp edges and tips are sometimes used with this device. Please store the device and all of its accessories and aids out of the reach of children. There is a risk of INJURY.

This device contains electronic components that are powered by either a mains connection or batteries. Never leave a child unsupervised with this device. The device should only be used as per these instructions otherwise there is a serious RISK of ELECTRICAL SHOCK.

Children should only use this device under supervision. Keep packaging materials (plastic bags, rubber bands, etc.) away from children. There is a risk of SUFFOCATION.

The chemicals and liquids provided should be kept out of reach of children. Do not drink the chemicals! Hands should be washed thoroughly under running water after use. In case of accidental contact with the eyes or mouth rinse with water. Seek medical treatment for ailments arising from contact with the chemical substances and take the chemicals with you to the doctor.

FIRE-/ DANGER OF EXPLOSION!

 Do not expose the device to high temperatures. Use only the mains adapter supplied or those battery types recommended. Never short circuit the device or batteries or throw into a fire. Exposure to high temperatures or misuse of the device can lead to short circuits, fire or even explosion!

RISK of material damage

 Never take the device apart. Please consult your dealer if there are any defects. The dealer will contact our service centre and send the device in for repair if needed.

Do not subject the device to temperatures exceeding 60° C.

WARNING! Not for children under three years. Choking hazard – small parts. Contains functional sharp edges and points!

ATTENTION: Only suitable for children of at least 8 years of age. Instructions for parents or other responsible persons are enclosed and must be followed. Keep packaging as it contains important information.

Check the Power supply for damage to the cord, plug, enclosure or other parts on a regular ba-

sis before use. Do not use until the damage has been repaired.

The toy is not intended for children under three years.

The toy must only be used with the recommended power supply

The power supply is not a toy.

The toy is not to be connected to more than the recommended number of power supplies.

Power supply Specification:

Rated voltage: 5V 

Input: 100-240V ~50/60Hz

Output: 5V 1000mA(DC)

TIPS on cleaning

Remove the device from its energy source before cleaning (remove plug from socket / remove batteries).

Clean the exterior of device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components.

Toys liable to be cleaned with liquid are to be disconnected from the power supply before cleaning

Clean the lens (objective and eyepiece) only with the cloth supplied or some other soft lint-free cloth (e.g. micro-fibre). Do not use excessive pressure - this may scratch the lens.

Dampen the cleaning cloth with a spectacle cleaning fluid and use it on very dirty lenses.

Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging. Batteries should be removed from the device if it is not going to be used for a long period of time.

DISPOSAL

Dispose of the packaging material/s as legally required. Consult the local authority on the matter if necessary.

Do not dispose of electrical equipment in your ordinary refuse. The European guideline 2002/96/EU on Electronic and Electrical Equipment Waste and relevant laws applying to it require such used equipment to be separately collected and recycled in an environment-friendly manner. Empty batteries and accumulators must be disposed of separately. Information on disposing of all such equipment made after 01 June 2006 can be obtained from your local authority.

EC Declaration of Conformity

Bresser GmbH has issued a „Declaration of Conformity“ in accordance with applicable guidelines and corresponding standards. The full text of the EU declaration of conformity is available at the following internet address:

www.bresser.de/download/8855000/CE/8855000_CE.pdf

Here are the parts of your microscope (Fig. 1-6):

- 1 10x WF Eyepiece
- 2 16x WF Eyepiece
- 3 Barlow Lens
- 4 Microocular
- 5 Eyepiece supports
- 6 Microscope Head
- 7 Set Screw
- 8 Objective Nosepiece
- 9 Objective
- 10 Clips
- 11 Microscope Stage
- 12 LED Illumination (transmitted light)
- 13 Microscope Base
- 14 Selection Knob for Illumination
- 15 Power supply
- 16 Focus knob
- 17 Color Filter
- 18 LED Illumination (reflected light)
- 19 Photomizer SE Software
- 20 5 Slides, 10 Cover Sips and 5 Prepared Specimens plastic box
- 21 Specimens:
 - a) Yeast
 - b) "Gum Media" Glue
 - c) Sea Salt
 - d) Shrimp Eggs
- 22 Transformer with Plug and Power Cable
- 23 Specimen Slicer
- 24 Shrimp Hatchery

- 25 Tweezers
- 26 Pipette
- 27 Carrying Case

How do I use my microscope?

Before you assemble your microscope, make sure that the table, desk or whatever surface that you want to place it on is stable, and does not wobble.

Apart from that, you will need a power source for your microscope's light (220V-330V).

How do I operate the electric LED illumination?

For the illuminator, you need the transformer with the plug and power cable (Fig. 6, 22). First, place the small plug on the power cable into the power connection on the back of the microscope (Fig. 1, 15). Then, insert the large plug into a power outlet.

There are two lights on the microscope. They do not use light bulbs, but rather light-emitting diodes (LED). The first lamp shines onto the specimen from below and the second from above. (The thing that you want to observe with the microscope is called the object or specimen, by the way.) You can use each lamp on its own, or both of them together. There is a selection knob for this (Fig. 1, 14). It has three numbers: I, II and III.

If you select the ...

- I, the light only comes from below (transmitted light).
- II, the light only comes from above (reflected light).
- III, both lamps shine light on the specimen.

For transparent objects (transmitted-light objects), number I is best. In order to observe firm, non-transparent objects (direct-light objects), select number II. For semi-transparent objects, it is best to select number III.

It is not recommended to use number III for transmitted-light objects on slides, since the light may cause reflections on the surface of the slide, which will disturb your observation.

When do I use the color filters?

The color filters (Fig. 1, 17) are located below the microscope stage (Fig. 1, 11). They help you when you are observing very bright or clear specimens. Here, you can choose from various colors. This helps you better recognize the components of colorless or transparent objects (e.g. grains of starch, protozoa).

How do I adjust my microscope correctly?

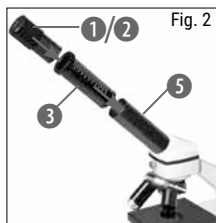
First, loosen the screw (Fig. 1, 7) and turn the microscope head (Fig. 1, 6) into a comfortable viewing position.

Each observation starts with the lowest magnification.

Adjust the microscope stage (Fig. 1, 11) so that it goes all the way down to the lowest position. Then, turn the objective nosepiece (Fig. 1, 8) until it clicks into place at the lowest magnification (objective 4x).

Note:

Before you change the objective setting, always move the microscope stage (Fig. 1, 11) to its lowest position. This way, you can avoid causing any damage!

Now, insert the 10x eyepiece (Fig. 1, 1) into the Barlow lens (Fig. 1, 3). Make sure that the Barlow lens is placed all the way into the eyepiece supports (Fig. 1, 5) and is not pulled out (Fig. 2).

How do I observe the specimen?

After you have assembled the microscope with

the adequate illumination and adjusted it correctly, the following basic rules are to be observed:

Start with a simple observation at the lowest magnification. This way, it is easier to position the object in the middle (centering) and make the image sharp (focusing).

The higher the magnification, the more light you will require for a good image quality.

Now place the prepared specimen (Fig. 6, 20) directly under the objective on the microscope stage (Fig. 3). The object should be located directly over the illumination (Fig. 1, 12).

In the next step, take a look through the eyepiece (Fig. 1, 1) and carefully turn the focus knob (Fig. 1, 16) until the image appears clear and sharp.

Now you can select a higher magnification by slowly removing the Barlow lens (Fig. 1, 3) from the eyepiece support (Fig. 1, 5). When the Barlow lens is almost completely pulled out, the magnification can be increased to almost double.

If you would like an even higher level of mag-

nification, insert the 16x eyepiece (Fig. 1, 2) and turn the objective nosepiece (Fig. 1, 8) to a higher setting (10x or 40x).

Important tip:

The highest magnification is not always the best for every specimen!

Note:

Each time the magnification changes (eyepiece or objective change, pulling out the Barlow lens), the image sharpness must be readjusted with the focus knob (Fig. 1, 16). When doing this, make sure to be careful. If you move the microscope stage too quickly, the objective and the slide could come into contact and become damaged!

Which light for which specimen?

With this unit, a reflected light and transmitted light microscope, you can observe transparent, semi-transparent as well as non-transparent objects.

The image of the given object of observation is "transported" through the light. As a result, only the correct light will allow you to see something!

If you are observing non-transparent (opaque) objects (e.g. small animals, plant components, stones, coins, etc.) with this microscope, the light falls on the object that is being observed.

From there, the light is reflected back and passes through the objective and eyepiece (where it gets magnified) into the eye. This is reflected light microscopy.

For transparent objections (e.g. protozoa), on the other hand, the light shines from below, through the opening in the microscope stage and then through the object.

The light travels further through the objective and eyepiece, where it is also magnified, and finally goes into the eye. This is transmitted-light microscopy.

Many microorganisms in water, many plant components and the smallest animal parts are already transparent in nature. Others have to be prepared. We may make them transparent through a treatment or penetration with the right materials (media), or by taking the thinnest slices from them (using our hand or a specimen slicer), and then examine them. You can read more about this in the following sections.

How do I make thin specimen slices?

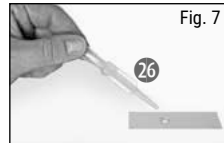
Only do this with the supervision of your parents or another adult.

As I already pointed out, the thinnest slices possible are taken from an object. In order to get the best results, we need some wax or paraffin. It is best if you get a candle. Place the

wax in a pot and heat it carefully over a low burner. Now, dip the object in the liquid wax a few times. Then, let the wax get hard. Using the specimen slicer (Fig. 6, 23) or a knife/scalpel, cut the smallest slices from the object that is covered with wax. These slices are to be laid on a slide and covered with a cover slip.

How do I make my own specimens?

Take the object that you want to observe and place it on a glass slide (Fig. 6, 20). Then, add a few drops of distilled water on the object (Fig. 7) using a pipette (Fig. 7, 26). Now, place a cover slip vertically at the edge of the drop of water, so that the water runs along the edge of the cover slip. Then, slowly lower the cover slip over the water drops (Fig. 8).

Note:

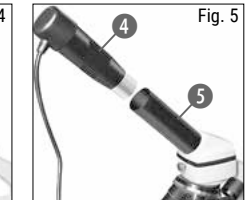
The included glue "gum media" (Fig. 5, 21b) is used to make permanent prepared specimens. Use this in place of the distilled water. If you want to keep the object in place on the slide permanently, use the gum media.

How do I install the MicOcular?

Note:

The MicOcular only works with the included Barlow lens! The magnification setting is changed when the MicOcular is used, and it must be readjusted with the focus knob.

First, remove the Barlow lens (Fig. 4, 3) with the current eyepiece from the eyepiece support (Fig. 4) and place the MicOcular (Fig. 5, 4) in the eyepiece (Fig. 5).

How do I install the software?

- 1 Insert the CD-ROM supplied in your CD or DVD drive. The installation menu opens automatically. If it doesn't, please switch to Windows Explorer and select your CD drive. Start the file "setup.exe" by double-clicking it.
2. A choice of languages is shown. Choose your preference. Confirm it by clicking „OK“. At the „Welcome“ screen, click „Next“>“.
3. In the next screen you will be asked for the „Destination Folder“. Confirm your input with

„Next>“.

- The setup status screen shows. This advises current operation status with a progress bar. This may take a few minutes.
- The „Completed the Photomizer Setup Wizard“ screen shows. Click here on „Finish“.

MikrOkular driver installation

During Photomizer SE software installation the correct driver for your operating system will be installed automatically. No manual input is needed.

In rare cases the device may not be recognised by your computer. As a rule you need then only install the driver manually from the CD. If this fails please refer to the troubleshooting chapter that follows.

Troubleshooting: the device isn't recognised by the computer (USB)

- This occurs in rare cases with some OEM Windows Vista or XP versions. The device is not defective. Updating the notebook or PC USB driver usually suffices.
- If not, proceed as below. (The maker cannot accept any liability whatsoever in these cases¹).

Delete ghost devices. These are devices not actually currently connected to your computer. The reason for this is that Windows generates a system settings entry for each USB device (e.g. USB Stick) on each port. It then seeks that/

those device/s each time it's started. This slows down USB device recognition and Windows booting and causes non-recognition of USB ports.

Cleaning up the system environment and deleting devices that are no longer in use helps here. To do so right click WORKPLACE (COMPUTER in Vista) and open the PROPERTIES menu. Then click the ADVANCED tab as in illustration 6 (ADVANCED SYSTEM SETTINGS in Vista) and then ENVIRONMENT VARIABLES. Choose the NEW option in ENVIRONMENT VARIABLES as in illustration 7. Enter the following under VARIABLE NAME as in illustration 8.

devmgr_show_nonpresent_devices

Under VARIABLE VALUE enter „1“.
Confirm your input with OK and then reboot the computer.

Change to the device manager. In the VIEW menu activate the SHOW DEVICES OPTION. The ghost devices previously not shown will now be displayed in light grey. Check the various categories including USB, memory size, etc. Only delete entries for devices you no longer use.¹

¹Source: Microsoft Knowledge Base
<http://support.microsoft.com/kb/315539/de>

*Not included

**System requirements: Intel X86 processor with at least 2.0 GHz speed, 200 MB hard disc space, 512 MB RAM memory, Windows XP or Vista (32-bit) and a free USB port (not via a hub if possible).

Using the MikrOkular

Preparation

- Slide a specimen under your microscope and focus on it.
- Remove the eyepiece and Barlow lens from the eyepiece support and the dust cap from your MikrOkular and install same in the eyepiece supports instead of the Barlow lens.
- Start your PC if you haven't yet and connect your MikrOkular to the USB port of your computer.

Showing and saving MikrOkular images on your PC.

- Start the Photomizer SE Software.
- Click „Open camera“
- If you have more than one device connected up select the one you want from the choice offered. Click „SoC PC-Camera“ or „MikrOkular“. Omit this step if you only have a single device connected up.
- The camera image should now be visible on your screen. Focus the microscope image.
- Click „Capture“ to record an image. It will then be shown on the right.
- Click on it to select it and then click „Transfer image“.
- This takes you to the Photomizer SE software.
- File - Save as

The Photomizer SE software

If you need help in the „Photomizer SE“ programme click „?“ and then „Open help“. If you need further assistance please visit the maker's home page at www.photomizer.net

Troubleshooting

Error	Solution
No recognizable image	<ul style="list-style-type: none">• Turn on light• Readjust focus
Image flickers (while observing with MicrOcular) = Monitor refresh rate not adequate	<ul style="list-style-type: none">• If necessary, increase resolution of graphics card

Make sure your microscope has a long service life.

Clean the lens (objective and eyepiece) only with the cloth supplied or some other soft lint-free cloth (e.g. microfibre). Do not press hard as this might scratch the lens.

Ask your parents to help if your microscope is really very dirty. The cleaning cloth should be moistened with cleaning fluid and the lens wiped clean using little pressure.

Make sure your microscope is always protected against dust and dirt. After use leave it in a warm room to dry off. Then install the dust caps and keep it in the case provided.

Warranty & Service

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at www.bresser.de/warranty_terms.

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